



# Purity Test of Epirubicin Hydrochloride with Reference to the Japanese Pharmacopoeia

Epirubicin Hydrochloride is a drug used in the treatment of various cancers. It is particularly important in the treatment of breast cancer.

A purity test was made for acetone, methanol, ethanol, 1-propanol, 1,4-dioxane (internal standard).

The standard was diluted with N,N-dimethylformamide. The resolution of acetone and 1,4-dioxane (internal standard) and the relative standard deviation of acetone, methanol, ethanol, and 1-propanol were calculated; excellent results were obtained.

## Example: Measurement of standard

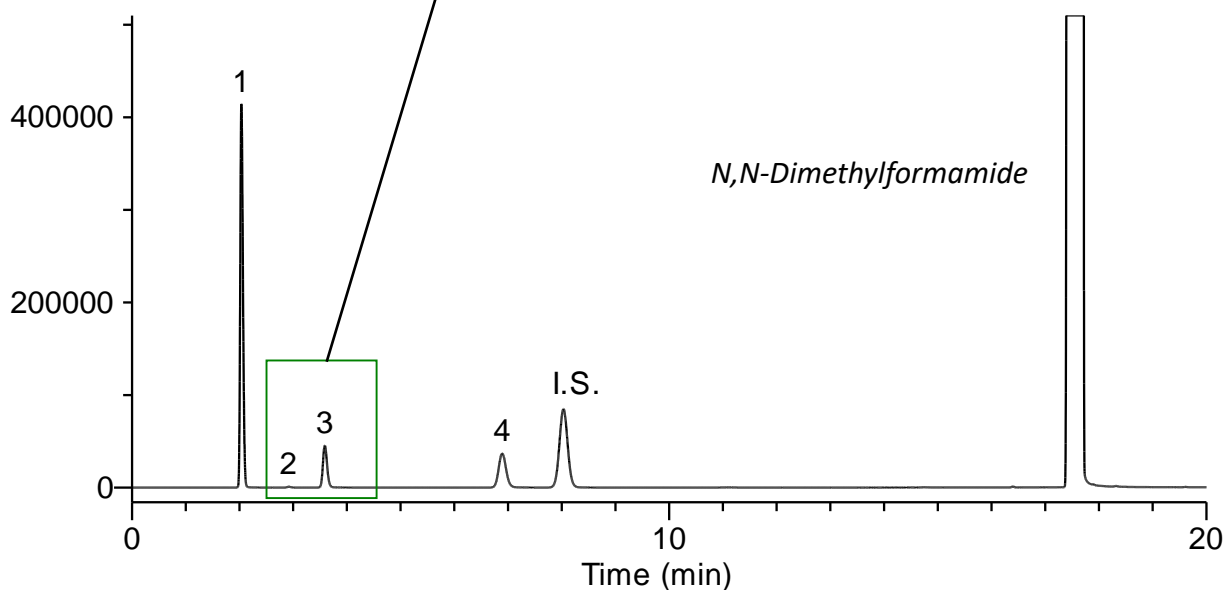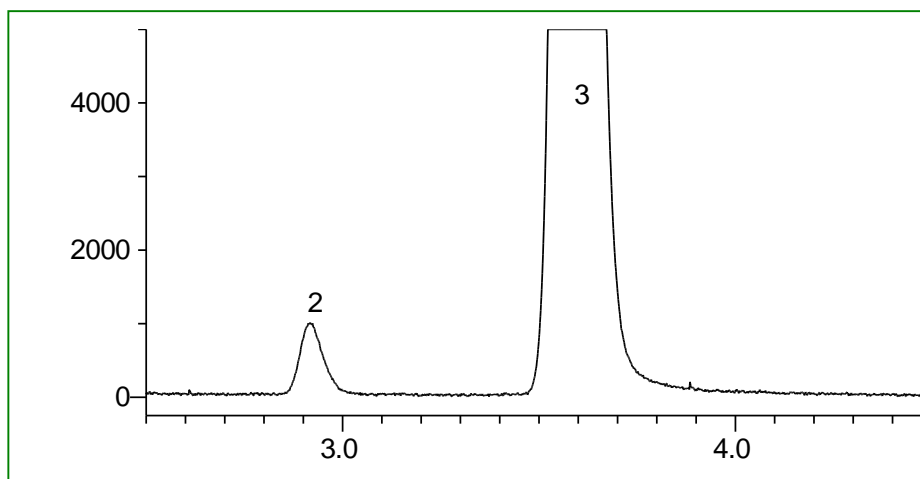

Analyte :

1. Acetone	1.25 $\mu\text{L}/\text{mL}$	4. 1-Propanol	0.32 $\mu\text{L}/\text{mL}$
2. Methanol	0.0068 $\mu\text{L}/\text{mL}$	I.S. 1,4-Dioxane	1.00 $\mu\text{L}/\text{mL}$
3. Ethanol	0.30 $\mu\text{L}/\text{mL}$		

\* The resolution between Acetone and 1,4-Dioxane is 32.1.

## Conditions

<b>System</b>	: GC - FID
<b>Column</b>	: InertCap WAX 0.53 mm I.D. × 30 m df = 1.00 μm
<b>Col. Temp.</b>	: 40 °C (11 min hold) - 10 °C/min - 90 °C - 50 °C/min - 130 °C (30 min hold)
<b>Carrier Gas</b>	: He 25 kPa
<b>Injection</b>	: Split 1:15 120 °C
<b>Detection</b>	: FID Range 10 <sup>0</sup> 150 °C
<b>Sample Size</b>	: Analyte in N,N-Dimethylformamide 1.0 μL

## Relative standard deviation

### Area repeatability (n = 6)

	Acetone	Methanol	Ethanol	1-Propanol	1,4-Dioxane (I.S.)
1 st	1398456	4260	241586	351661	899222
2 nd	1396842	4272	241729	352627	901190
3 rd	1394689	4300	240834	351744	897514
4 th	1393229	4209	240332	350138	895809
5 th	1385342	4111	239718	349265	891042
6 th	1383033	4085	239039	349311	890195
Ave.	1391932	4206	240540	350791	895829
SD	6302	89	1055	1413	4422
RSD (%)	0.5	2.1	0.4	0.4	0.5

GL Sciences disclaims any and all responsibility for any injury or damage which may be caused by this data directly or indirectly. We reserve the right to amend this information or data at any time and without any prior announcement.

### GL Sciences, Inc. Japan

22-1 Nishishinjuku 6-Chome  
Shinjuku-ku, Tokyo,  
163-1130, Japan  
Phone: +81-3-5323-6620  
Fax: +81-3-5323-6621  
Email: [world@glsciences.com](mailto:world@glsciences.com)  
Web: [www.glsciences.com](http://www.glsciences.com)

### GL Sciences B.V.

De Sleutel 9  
5652 AS Eindhoven  
The Netherlands  
Phone: +31 (0)40 254 95 31  
Email: [info@glsciences.eu](mailto:info@glsciences.eu)  
Web: [www.glsciences.eu](http://www.glsciences.eu)

### GL Sciences, Inc. USA

4733 Torrance Blvd. Suite 255  
Torrance, CA 90503  
Phone: 310-265-4424  
Fax: 310-265-4425  
Email: [info@glsciencesinc.com](mailto:info@glsciencesinc.com)  
Web: [www.glsciencesinc.com](http://www.glsciencesinc.com)

### GL Sciences (ShangHai) Ltd.

Tower B, Room 2003,  
Far East International Plaza,  
NO,317 Xianxia Road,  
Changning District.  
Shanghai, China P.C. 200032  
Phone: +86 (0)21-6278-2272  
Email: [contact@glsciences.com.cn](mailto:contact@glsciences.com.cn)  
Web: [www.glsciences.com.cn](http://www.glsciences.com.cn)

### International Distributors

Visit our Website at [www.glsciences.com/distributors](http://www.glsciences.com/distributors)