# Analysis of cyclosporin A in serum using MonoSpin C18 and Ph.

LT203 GL Sciences Inc.

Cyclosporin A (Cyclosporin A) is a type of cyclic polypeptide antibiotic and is used as a medical drug. It is used to suppress the immune system of organ transplant patients, thereby reducing the risk of organ transplant rejection. However, due to the narrow effective therapeutic window and the potential for toxicity, it is necessary to monitor blood drug concentrations. In this study, cyclosporin A was added to human serum, extracted and purified using spin columns Monospin C18 and Ph, and analyzed by LC-MS/MS. (T.Kunieda)

#### 1. Pretreatment with MonoSpin

#### ♦ What is MonoSpin?

The MonoSpin series is a spin column using a silica monolith with uniform continuous pores. Silica monolith with high porosity is used as a carrier, so liquid can be passed through only by centrifugation. Therefore, it is possible to purify and concentrate the sample with a short and simple operation. In addition, the bed volume is small and the liquid is easily drained, so it is also suitable for small sample volumes. 1

MonoSpin



Monolithic disc

how to use  $\rightarrow$  <u>https://youtu.be/YDhZ1Dtwq3k</u>

application  $\rightarrow$  <u>https://www.gls.co.jp/technique/lifescience/monospin/application/index.html</u>



#### sample solution

uman serum: 0.1% formic acid: cyclosporin (50% acetonitrile containing 0.1% formic acid) = 200:190:10µL, centrifuged at 10,000 x g for 1 min, and the supernatant was used as the sample solution. \* By removing fine particles by centrifugation in advance, the flow of the spin column will be improved.

		Rec (%)	RSD(%)
0.5 ng/mL	C18	94.8	2.4
	Ph	103.8	10.0
5 ng/mL	C18	98.6	12.1
	Ph	115.6	14.8

#### Spike recovery test results

(0.5,5 ng/mL in sample, n=3 x 1 days)

Cyclosporin A



1.5e5

1.4e5 1.3e5 1.2e5 1.1e5

7 0ea

5.0e4

4.0e4

3.0e4

2.0e4

0.0

0.0

Chromatogram example of standard solution



	Time (min)	A (VOI%)	B (VOI%)
	0.0	10	90
	7.0	90	10
	10.0	90	10
	10.1	10	90
	13.0	10	90
:	0.4 mL/min		

0 0 05 10 1.5 20 2.5 3.0 3.5 4.0 4.5 5.0 5.5 6.0 6.5 7.0 7.5 8.0 8.5 9.0 9.5 10.0 10.5 11.0 11.5

## Examination of LC column and analysis conditions

#### ■ InertSustain AQ-C18 (1.9 µm 100 x 2.1 mm I.D.)

InertSustainSwift C8 (3 μm 100 x 3.0 mm I.D.)

5.0

Since the target component is highly hydrophobic, an initial study was conducted with an initial concentration of acetonitrile of 50% and an oven temperature of 40°  $\,$  C.

<sup>5.0</sup> Time, min

Time, min

10.0

10.0

Flow Rate

Since the retention was strong with the ODS column and the peak was broadened, we investigated with the C8 column.



Since the peak width was wide even with the C8 column, we next investigated with the phenyl column.





The phenyl column gave narrower peak widths, and

changing the eluent gradient resulted in sharper peaks. The peak shape was further improved by increasing the column oven temperature to 50°C.

#### **3.Related product**

#### **InertSustain Phenyl**

Phenyl-bonded columns give a different separation pattern than C18 columns due to  $\pi$ electron interactions. It is effective as a column of choice when separation is insufficient with a C18 column. In particular, InertSustain Phenyl has a phenyl group directly bonded to silica gel, so it has the characteristic of recognizing the difference in the electronic state of aromatic compounds to a greater extent than a general phenyl column (alkylphenyl groupbonded columns). In addition, since the phenyl group is chemically modified at high density, it has excellent stereoselectivity and exhibits unprecedented separation performance.



Column	Size	Cat.No.
InertSustain Phenyl	3µm, 50 x 2.1 mm I.D.	5020-16433

### MonoSpin(Solid-phase extraction column for minute samples)

MonoSpin is a pretreatment spin column that uses a silica monolith with uniform continuous pores. All operations can be performed with a short centrifugation process.

The S type is ideal for samples up to 800  $\mu$ L, the L type is ideal for larger 1-8 mL samples, and the 96-well plate type is available for multi-sample processing.



#### MonoSpin C18

A column with octadecyl groups and reversed-phase partitioning interaction. Ideal for extracting drugs from biological samples and desalting/concentrating peptide samples.

Product name	Quantity	Cat.No.
MonoSpin C18	50	5010-21700



MonoSpin S type

#### **MonoSpin Ph**

Reversed-phase mode type with chemically modified phenyl group. It is possible to perform pretreatment with higher selectivity by utilizing the weaker hydrophobic interaction than C18. Suitable for recovering hydrophobic drugs from biological samples.

Product name	Quantity	Cat.No.
MonoSpin Ph	50	5010-21733

GL Sciences disclaims any and all responsibility for any injury or damage which may be caused by this data directly or indirectly. We reserve the right to amend this information or data at any time and without any prior announcement.

GL Sciences Inc. Japan 22-1 Nishishinjuku 6-chome Shinjuku-ku, Tokyo 163-1130, Japan

 Phone: +81-3-5323-6620

 Fax:
 +81-3-5323-6621

 Email:
 world@gls.co.jp

 Web:
 www.glsciences.com



International Distributors Visit our Website at www.glsciences.com/distributors

GL Sciences Inc. USA 4733 Torrance Blvd. Suite 255 Torrance, CA 90503 USA

 Phone:
 +1-310-265-4424

 Fax:
 +1-310-265-4425

 Email:
 info@glsciencesinc.com

 Web:
 www.glsciencesinc.com

**GL Sciences B.V.** Dillenburgstraat 7C 5652AM, Eindhoven The Netherlands

Phone: +31-40-254-9531 Email: info@glsciences.eu Web: www.glsciences.eu

#### **GL Sciences (Shanghai) Limited**

Tower B, Room 2003 Far East International Plaza No.317 Xianxia Road, Changning District Shanghai, China 200051

Phone: +86-21-62782272 Email: contact@glsciences.com.cn Web: www.glsciences.com.cn