


Analysis of Hydrocarbons C5-C10

Specification

One GC Channel

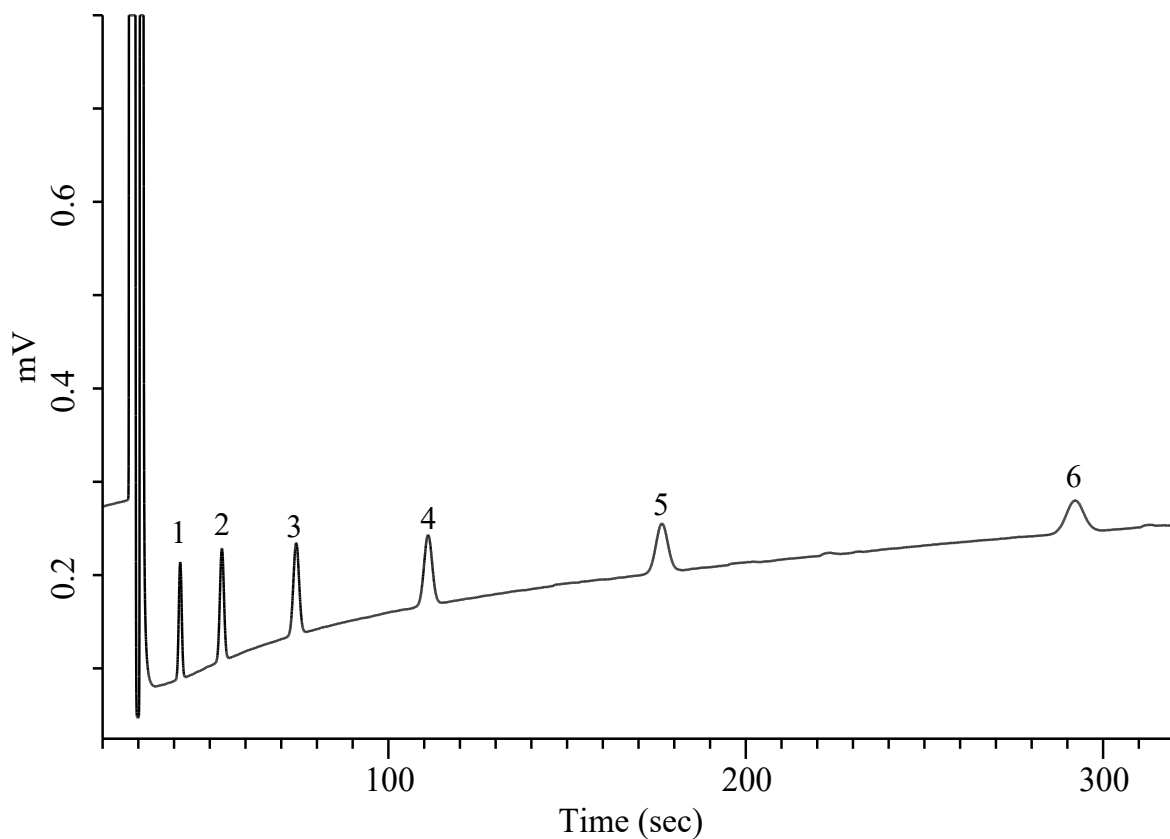

Conditions

System : Micro GC
Column : CP-Volamine 15 m
Column Temp. : 120 °C
Carrier Gas : He 190 kPa
Inject Time : 40 msec

Analyte:

1. <i>n</i> -Pentane (C ₅ H ₁₂)	50 ppm (v/v)
2. <i>n</i> -Hexane (C ₆ H ₁₄)	50 ppm (v/v)
3. <i>n</i> -Heptane (C ₇ H ₁₆)	50 ppm (v/v)
4. <i>n</i> -Octane (C ₈ H ₁₈)	50 ppm (v/v)
5. <i>n</i> -Nonane (C ₉ H ₂₀)	50 ppm (v/v)
6. <i>n</i> -Decane (C ₁₀ H ₂₂)	50 ppm (v/v)

Balance Gas Air